

Femtomole Peptide Analysis by PicoFrit® Nanobore LC/MS at Low Pressures

Introduction

The inherent chemical specificity and sensitivity of electrospray ionization mass spectrometry (ESI-MS) has led to the development of integrated nanoscale liquid chromatography (nLC) ESI systems.^{1,2} In this approach, an appropriate ESI emitter is fabricated directly on the nLC column outlet. Method development has been severely limited by the difficult fabrication of suitable integrated nLC-ESI columns. Furthermore, instrumentation necessary for the generation of suitable sample injection and subsequent chromatography has been specialized and complex. The combination of a tapered, fritted fused-silica needle packed with a highporosity reverse-phase media eliminates these difficulties. This device, the PicoFrit®, provides purification, concentration, and separation at low column pressures. Low-pressure operation eliminates the need for specialized HPLC hardware, provides for short nLC run times, and allows direct integration with common syringe pumps and/or auto-samplers as shown in Figure 1.

Column Fabrication

PicoFrit® columns were fabricated from fused-silica tubing with a 30 µm ID tip, an integral high-porosity frit, and multi-layer conductive coating (PF360-75-30-CE) as shown in Figure 2. Columns were syringe-packed with 10 µm POROS®, R2 phase media as follows: Approximately 200 ml of freshly ultrasonicated POROS slurry in MeOH (5 mg/ml) was drawn into a 500-ml gas-tight Luer-lock syringe (Hamilton Company). The distal end of a 50-cm ESI column was inserted into the barrel of the syringe using Luer/fused-silica adapter components from Upchurch Scientific®. Columns were packed by hand pressure alone; packing progress was monitored by light microscopy. When the desired length (about 5 cm) was reached, the column was rinsed with 50 ml of MeOH. Columns were dried for long-term storage with dry nitrogen at 500 psi for 15 minutes. Prior to use, columns were re-hydrated with MeOH and equilibrated with 1% acetic acid.

Advantages of PicoFrit® Combined Column/Emitter

- Provides routine high-sensitivity, low-fmol limit of detection, for peptides on an ion trap in MS/MS mode
- Dirty and/or dilute peptide mixtures now easily run in ESI mode
- Zero post-column effects (e.g. sample loss, resolution loss)
- Capable of fast sample turn-around (< 5 min/run) for high throughput
- Packing the column in the tip eliminates problems with clogged tips
- Operable at low (syringe pump) pressures, eliminating the need for specialized hardware

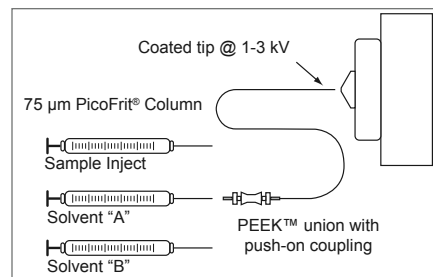


FIGURE 1 Linking a hand inject, rinse or syringe pump and a PicoFrit® on a Thermo Finnigan LCQ™

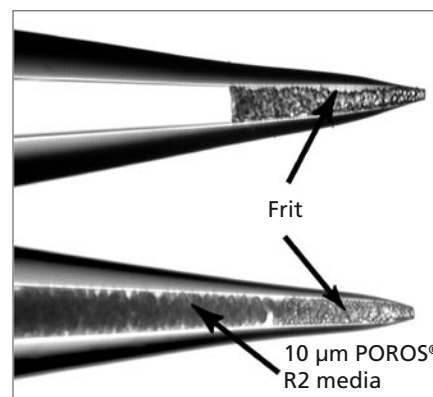


FIGURE 2 Unpacked (top) and pre-packed (bottom) PicoFrit® columns

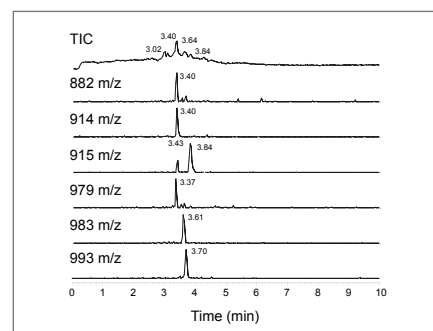


FIGURE 3 75 m x 47 mm (200 nL) POROS® R2, 30 m PicoFrit® @ 1 L/min 30:70:1 (ACN:H₂O:HOAc) (10 L inj. 79 - 165 fmol/peptide)

nLC/ESI-Mass Spectrometry

LC-ESI columns were mounted on a LCQ™ ion-trap mass spectrometer (Thermo Finnigan Inc.) using an inline tip adapter, model ADPT-TLC (New Objective, Inc.) shown in Figure 4. High-voltage (1-3 kV) contact was established with the coated end of the needle using a spring contact mechanism. The distal end of the column was attached to a modified PEEK™ union (Upchurch Scientific). The other end of the union was fitted with a tubing sleeve that provided for rapid push-on coupling to a variety of stainless-steel needles on gas-tight syringes (SGE Inc.). On-column sample injection, washing, and elution were performed by manually swapping syringes. Injection was performed by hand, while solvent flow for washing and elution (0.2 to 1 ml/min) utilized the LCQ syringe pump. MS/MS data was acquired in a data-dependent manner. A test sample consisting of a mixture of six synthetic class I kb murine peptides (7 to 8 bases long, covering a mass range of 881 to 992 Da) was prepared in 1% acetic acid, final concentrations ranging from 7.9 to 16.5 fmol/ml per peptide.



FIGURE 4 ADPT-TLC on the Thermo Finnigan LCQ™

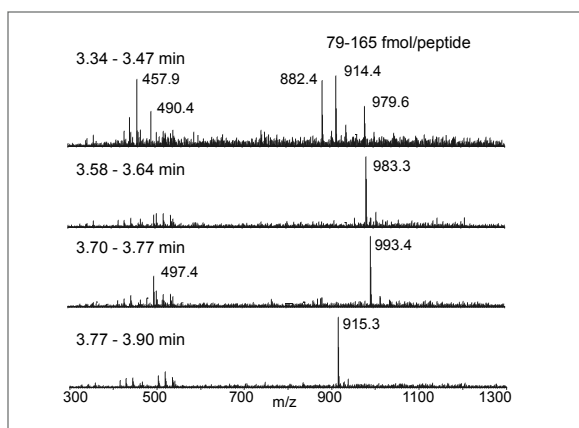


FIGURE 5 MS Results

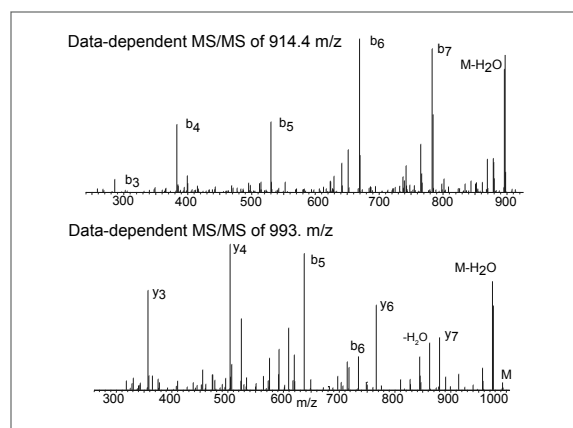


FIGURE 6 MS/MS Results

MS/MS Results Acknowledgments

For the LCQ™ data: Andrew J. Tomlinson and Roman M. Chicz, Pangaea Pharmaceuticals, Lexington, MA. For supplying the synthetic peptide samples: Steve Jameson and Chad Kieper, Center for Immunology, University of Minnesota Medical School.

References

- 1) M. R. Emmett and R. M. Caprioli. "Micro-Electrospray Mass Spectrometry: Ultra-High-Sensitivity Analysis of Peptides and Proteins," *J. Am. Soc. Mass Spectrom.* 1994, 5, 605-613
- 2) M. T. Davis, D. C. Stahl, S. A. Hefta, T. D. Lee. "A Microscale Electrospray Interface for On-Line Capillary Liquid Chromatography/Tandem Mass Spectrometry of Complex Peptide Mixtures." *Anal. Chem.* 1995, 67, 4549-4556

The information contained in this circular is believed reliable and accurate; however, nothing set forth herein constitutes a warranty or representation of any kind or nature. Caution: Particular end-user applications for these products may be restricted by existing patents. Complying with any such patent is the sole responsibility of the user. PicoFrit is a registered trademark of New Objective, Inc. POROS is a registered trademark of Applied Biosystems. LCQ is a trademark of the Thermo Electron Corporation. PEEK is a trademark of Victrex plc. New Objective, Inc. reserves the right to change specifications of its products without notice. PicoFrit columns are manufactured under U.S. Patents 5,997,746 and 6,190,559 and sold under license of U.S. Patent 5,572,023. © 2004 New Objective, Inc. All rights reserved.